



9520 Series Pulse Generator  
Operating Manual

QUANTUM COMPOSERS, INC  
PO Box 4248  
Bozeman, MT 59772  
(406)582-0227 phone  
(406)582-0237 fax  
[www.quantumcomposers.com](http://www.quantumcomposers.com)

# Contents

<b>1 Introduction .....</b>	<b>1</b>
Technical Support .....	2
Warranty .....	2
Package Contents .....	2
Safety Issues .....	3
<b>2 Front Panel Overview .....</b>	<b>5</b>
<b>9520 Front Panels .....</b>	<b>6</b>
Display Layout and Indicators .....	6
<b>Description of Front-Panel Area .....</b>	<b>6</b>
Keypads .....	6
Rotary Adjustment Knob .....	7
Second Level Menus .....	7
<b>3 Pulse Concepts &amp; Pulse Generator Operation .....</b>	<b>9</b>
Counter Architecture Overview .....	10
System Timer Functions .....	10
Channel Timer Functions .....	11
Digital Output Multiplexer .....	11
Dependent & Independent Timing Events .....	11
<b>Navigating the 9520 Front Panel .....</b>	<b>11</b>
Selecting Menus .....	11
Selecting Menu Items .....	12
Numeric Input Mode .....	12
Entering Non-Numeric Parameters .....	12
Alphanumeric Input Mode .....	12
Enabling System Output .....	13
Enable/Disable Channel Output .....	13
Rearming the Channel Timers .....	13
Setting Pulse Timing Parameters .....	13
Setting Pulse Output Parameters .....	13
Using the Output Multiplexer .....	14
Setting System Internal Rate Parameters .....	14
<b>4 9520 Menus .....</b>	<b>15</b>
<b>9520 Menu Structure .....</b>	<b>16</b>
<b>System Mode Menu .....</b>	<b>18</b>
Setting System Mode of Operation .....	18
<b>Channel Menus .....</b>	<b>18</b>
Enabling Channel Output .....	19
Setting the Channel Timing Parameters .....	19
Setting Pulse Output Configuration .....	19
Setting Channel of Operation .....	19

Delaying the Start of Channel Output .....	19
Configuring the Channel Multiplexer .....	19
<b>Advanced Channel Menus .....</b>	<b>20</b>
Setting the Sync Source .....	20
Setting Channel Gate Control .....	20
<b>Gate Menu .....</b>	<b>20</b>
Enabling System Gate .....	21
<b>Trig Menu .....</b>	<b>21</b>
Enabling System Trigger .....	21
<b>Rate Menu .....</b>	<b>21</b>
Setting the Clock Source and Internal Rate .....	21
Setting the Output Reference .....	21
<b>System Menu .....</b>	<b>22</b>
Setting System Communication Parameters .....	22
Setting Keypad Parameters .....	22
Setting the Auto Start Mode .....	22
Setting the Display Decimal Mark .....	22
Setting the Display Brightness .....	22
<b>Store Menu .....</b>	<b>23</b>
Storing a Configuration .....	23
<b>Recall Menu .....</b>	<b>23</b>
Recalling System Configurations .....	23
<b>Counter Menu .....</b>	<b>24</b>
Using the Counter Function .....	24
<b>Information Menu .....</b>	<b>24</b>
<b>5 Operating the 9520 .....</b>	<b>25</b>
<b>Quick Start - Normal Internal Rate Generator Operation .....</b>	<b>26</b>
<b>Quick Start - Normal External Trigger Operation .....</b>	<b>26</b>
<b>System Timer Overview .....</b>	<b>27</b>
To Use Continuous Mode .....	27
To Use Single Shot Mode .....	27
To Use System Burst Mode Function .....	28
To Use System Duty Cycle Function .....	28
<b>Channel Timer Overview .....</b>	<b>28</b>
To Use Normal Mode Function .....	28
To Use Channel Single Shot Function .....	29
To Use Channel Burst Mode Function .....	29
To Use the Channel Duty Cycle Function .....	29
To Use the Channel Gating Function .....	30
<b>External Input Overview .....</b>	<b>30</b>
To Generate a Pulse on Every Trigger Input .....	30
To Generate a Burst of Pulses on Every Trigger Input .....	31
To Start a Continuous Stream of Pulses Using the External Trigger .....	31
To use the External Gate to Control the System .....	32
<b>6 Programming the 9520 .....</b>	<b>33</b>
<b>Personal Computer to Pulse Generator Communication .....</b>	<b>34</b>

RS232 Interface Overview .....	34
USB Interface Overview .....	34
GPIB Interface Overview .....	35
Ethernet Interface Overview .....	35
Programming Command Types and Format .....	36
Line Termination .....	36
IEEE 488.2 Common Command Format .....	36
SCPI Command Keywords .....	37
SCPI Command Format .....	37
SCPI Format .....	37
SCPI Keyword Separator .....	37
SCPI Optional Keywords .....	38
SCPI Specific and Implied Channel .....	38
SCPI Parameter Types .....	38
Error Codes .....	38
Programming Examples .....	39
9520 SCPI Command Summary .....	40
9520 SCPI Command Summary .....	41
9520 SCPI Command Summary .....	42
9520 SCPI Command Summary .....	43
IEEE 488.2 Common Commands .....	44
<b>Appendix .....</b>	<b>45</b>
Appendix A - 9520 Specifications .....	46
Appendix B - Safety Symbols .....	49
Appendix C - 35V Output Module .....	53
Appendix D - Dual Trigger Input .....	54
Appendix E - Impedance Matching Outputs .....	56
Appendix F - Com.....	57
Appendix G - Ethernet Connectivity .....	59
Appendix H - Increment Mode .....	61
Appendix I - SRS Command Set Compatibility .....	63
Appendix J - AT45 Volt Module .....	64



---

# Introduction

This manual is a reference designed to familiarize you with the Quantum Composers 9520 series pulse generator and is arranged so that you can easily find the information you're looking for. Generally, each topic has its own section and no section assumes that you've read anything else in the manual.

## Technical Support

For questions or comments about operating the 9520 -- our technical staff can be reached via one of the following methods:

- Phone - (406) 582-0227
- Fax - (406) 582-0237
- Internet - [www.quantumcomposers.com](http://www.quantumcomposers.com)

## Warranty

In addition to a 30-day money back guarantee, the 9520 has a two-year limited warranty from the date of delivery. This warranty covers defects in materials and workmanship. Quantum Composers will repair or replace any defective unit. Contact us for information on obtaining warranty service.

## Package Contents

The box you receive should contain the following:

- 9520 Pulse Generator
- AC Power Cord
- User's Manual on Disk

Contact Quantum Composers (406) 582-0227 if any parts are missing.

## Safety Issues

Normal use of test equipment presents a certain amount of danger from electrical shock because testing must be performed where exposed voltage is present.

An electrical shock causing 10 milliamps of current to pass through the heart will stop most human heartbeats. Voltage as low as 35 VDC or RMS AC should be considered dangerous and hazardous since it can produce a lethal current under certain conditions. Higher voltages pose an even greater threat because such voltage can easily produce a lethal current. Your normal work habits should include all accepted practices that will prevent contact with exposed high voltage, and steer current away from your heart in case of accidental contact with a high voltage. You will significantly reduce the risk factor if you know and observe the following safety precautions:

- If possible, familiarize yourself with the equipment being tested and the location of its high-voltage points. However, remember that high voltage may appear at unexpected points in defective equipment.
- Do not expose high voltage needlessly. Remove housing and covers only when necessary. Turn off equipment while making test connections in high-voltage circuits. Discharge high-voltage capacitors after shutting down power.
- When testing AC powered equipment, remember that AC line voltage is usually present on power input circuits, such as the on-off switch, fuses, power transformer, etc.
- Use an insulated floor material or a large, insulated floor mat to stand on, and an insulated work surface on which to place equipment. Make certain such surfaces are not damp or wet.
- Use the time-proven “one hand in the pocket” technique while handling an instrument probe. Be particularly careful to avoid contact with metal objects that could provide a good ground return path.
- Never work alone. Someone should always be nearby to render aid if necessary. Training in CPR first aid is highly recommended.



---

**Front Panel Overview**

---

---

# Front Panel Overview

## 9520 Front Panels

### Display Layout and Indicators

A 4 line x 20 character vacuum fluorescent display module displays parameters and status information. The status information is located in the upper-left corner of the display, between the two brackets. There are four enunciators:

- Vertical Arrow                      Indicates there are additional pages to the current menu.
- Blinking Light                      Indicates the unit is actively generating pulses, or armed and waiting for an external trigger.
- Musical Note                        Indicates the function key has been pressed.
- Question Mark                      In external oscillator operation, indicates the internal pll is not yet locked with the external clock signal.

The upper-right side of the display contains the title of the currently displayed menu. The rest of the display is used for system parameters. The display brightness may be adjusted, allowing the instrument to be used under various lighting conditions.

### Description of Front-Panel Area

#### Keypads

Three keypad areas provide fast access to various menus and easy editing of system parameters.

- Channel Keypad                      Provides one touch access to the menus for setting up the channel parameters. Pressing the appropriate letter will display the parameters for the corresponding channel.
- Arrow Keypad                        The up/down arrows are used to increment/decrement the current parameter (indicated by the blinking cursor). The position of the cursor controls the step size for each increment. The right/left arrow moves the cursor to different positions within the current parameter. The **NEXT** key selects the next parameter in the currently displayed menu.
- Numeric Keypad                      Allows numbers and alphanumeric values to be entered. When entering alphanumeric values, pressing a key will display the first letter shown on the key. The yellow **FUNC** key allows the keys to select the yellow functions.

Repeated key presses will toggle through all the letters, both upper and lower case, shown on the keycap. To enter two letters which appear on the same keycap, select the first character, then use the right arrow to shift to the next position and enter the next letter. When data entry is complete the **ENTER** key must be pressed.

### **Rotary Adjustment Knob**

An alternate to the Arrow Keypad, the Rotary Adjustment Knob may be used to adjust the current parameter. The step size is controlled by the position of the cursor, however turning the knob faster will increase the step size. Pushing the knob will perform functions similar to the **NEXT** key and switch to the next parameter in the currently displayed menu.

### **Second Level Menus (Function Key)**

The second level menus (indicated in yellow above certain keys) are accessed through the use of the yellow **FUNC** key. Pressing the **FUNC** key once and then pressing the desired menu key will display the specified second level menu. Pressing the **FUNC** key twice in succession will put the unit into “Function Lock” mode, where the second level menus can be accessed without repeatedly pressing the **FUNC** key. Pressing the **FUNC** key a third time will exit “Function Lock” mode.



---


**Pulse Concepts and  
Pulse Generator Operations**

---

---

# Pulse Concepts and Pulse Generator Operation

## Counter Architecture Overview

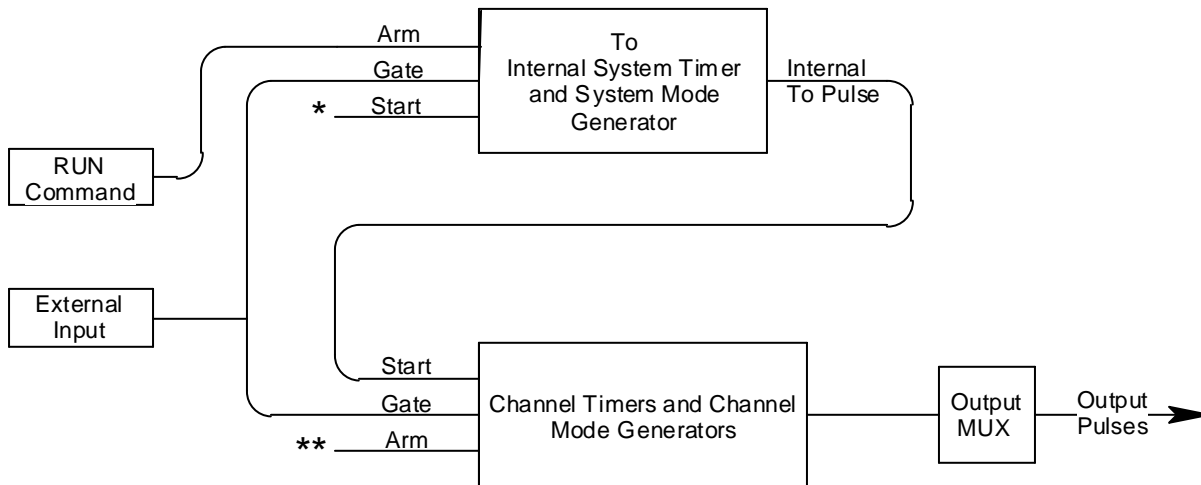

\*Start source is: RUN button in Internal Modes  
External input in External Trigger modes  
\*TRG command via Serial/GPIB access

\*\*Channels are armed by the RUN button. In single shot and burst modes channels may be rearmed by pressing the RUN button.

## System Timer Functions

The System Timer functions as a non-retriggerable, multi-vibrator pulse generator. This means that once started, depending on the mode, the timer will produce pulses continuously. Before pulses can be generated, the timer must be armed and then receive a start pulse. Arming the counter is done by pressing the **RUN/STOP** key. With external trigger disabled, the **RUN/STOP** key also generates the start command for the counter. With external trigger enabled, the external trigger provides the start pulse. In either case, once started, the counter operation is determined by the System Mode Generator. Standard modes include:

- Continuous                      Once started  $T_0$  pulses are generated continuously.
- Single Shot                      One  $T_0$  pulse is generated for each start command.
- Burst                              'n'  $T_0$  pulses are generated for each start command.
- Duty Cycle                      Once started  $T_0$  pulses cycle on and off continuously.

The  $T_0$  pulses are distributed to all of the start inputs of the Channel Timers and Mode Generators

## Channel Timer Functions

The Channel Timer functions as a non-retriggerable, delayed, one shot pulse generator. This means that the timer will only generate one delayed pulse for every start pulse received. Once the channel timer has started counting, additional start pulses will be ignored until the pulse has been completed (non-retriggerable). The start pulse for each channel is provided by the internal  $T_0$  pulse generated by the Internal System Timer. Whether or not a pulse is generated for each  $T_0$  pulse is determined by the Channel Mode Generator. Standard modes include:

- Normal A pulse is generated for each  $T_0$  pulse.
- Single Shot One pulse is generated at the first  $T_0$  pulse, after which output is inhibited.
- Burst A pulse is generated for each  $T_0$  pulse, 'n' times, after which output is inhibited.
- Duty Cycle 'n' pulses are generated for each  $T_0$  pulse after which the output is inhibited for 'm' times. The cycle is then repeated.

Different modes may be selected for each output, allowing a wide variety of output combinations. Each output may also be independently disabled or gated (using the external gate input).

## Digital Output Multiplexer

The output of the Channel Timers are routed to a set of multiplexers. This allows routing of any or all Channel Timers to any or all of the unit outputs. In the normal mode of operation, the output of the nth Channel Timer is routed to the Tnth output connector. As an example, if a double pulse is required on Channel A output, one can multiplex the Channel A timer with the Channel B timer adjusting each timer to provide the necessary pulses.

## Dependent & Independent Timing Events

The 9520 allows the user to control the relationship between the Channel Timers by setting the sync source for each timer. Independent events are all timed relative to the internal  $T_0$  start pulse. Dependent events may be linked together by setting the sync source to the controlling event. This allows the instrument to match the timed events and adjustments can be made in one event without detuning the timing between it and a dependent event.

## Navigating the 9520 Front Panel

### Selecting Menus

Parameters are grouped in menus, selectable using the function keys. To select the output channel parameters press the letter key corresponding to the desired channel. To select other menus, including the advanced channel menus, press the **FUNC** key and then the key corresponding to the desired function.

Menus may include a number of different pages. Each page containing up to four parameters. The status block in the upper-left corner of the display shows a vertical arrow if the current menu contains additional pages. To select the next page, press the channel button again or select the same menu pressing the **FUNC** key and the channel/menu key again.

### Selecting Menu Items

Within a menu, the blinking cursor indicates the current menu item for editing. The **NEXT** key or pressing the adjustment knob will select a different menu item.

### Numeric Input Mode

When the current item is numeric, the system enters the Numeric Input Mode. In this mode data may be edited in one of three ways. Using the arrow keypad, the Left and Right arrow keys are used to select a digit to edit. The selected digit blinks to identify itself as the active digit. The Up and Down arrow keys are then used to increment or decrement this digit. Alternately, after using the Left and Right arrow keys to select an active digit, the adjustment knob may be used to increment and decrement this digit. The adjustment knob features speed dependent resolution. Slow rotation will increment or decrement the active digit by one. As you increase the speed of rotation, the parameter will be 10 to 1000 times faster depending on the speed.

The last entry mode is using the numeric keypad. Enter the number, including decimal point using the numeric keypad. Complete the number using the **ENTER** key. Errors may be corrected using the backspace key or to start over press the clear key (**CLR**). Pressing the **CLR** key a second time will exit the numeric keypad mode and restore the original number.

### Entering Non-Numeric Parameters

When the current item is non-numeric, the Up and Down arrow keys are used to select among different options for the parameter. The adjustment knob may also be used to change the selection. If the item is an on-off toggle, the Up arrow (CW adjustment knob) enables the item and the Down arrow (CCW adjustment knob) disables the item.

### Alphanumeric Input Mode

When the current item is alphanumeric, the system enters the Alphanumeric Input Mode. In this mode, data is entered using the alphanumeric keypad. When entering alphanumeric values, pressing a key will display the first letter shown on the keypad. Repeated key presses will toggle through all the letters, both upper and lower case, shown on the key cap. To enter two letters which appear on the same key cap, select the first character, then use the right arrow to shift to the next position and enter the next letter. The Left and Right arrow keys may be used to position the cursor to edit any character. When data entry is complete, the **ENTER** key must be pressed. The keys contain the following characters:

- 1                    1 2 3 4 5 6 7 8 9 0
- 2                    A B C a b c 2

- 3 DEFdef3
- 4 GHIghi4
- 5 JKLjkl5
- 6 MNOmno6
- 7 PQRSpqrs7
- 8 TUVtuv8
- 9 WXYZwxyz9
- 0 0123456789
- . , # \$ % & ?
- - + \* / space

### Enabling System Output

The RUN/STOP key is used to arm the system. With external trigger disabled, the key will arm and start pulse output. With external trigger enabled, the key will arm the pulse generator. Pulse output then starts after the first valid trigger input. Pressing the RUN/STOP key a second time disables the pulse generator.

### Enable/Disable Channel Output

At the top of each channel menu page is a parameter to enable or disable the output of the channel. Each channel may be individually enabled or disabled. An illuminated channel key indicates that the channel is enabled.

### Rearming the Channel Timers

In the channel single shot mode and burst mode, the Channel Timers may be rearmed after completing the initial output by pressing the **FUNC** key followed by the **RUN/STOP** key. If there are channels currently running in normal mode, single shot and burst channels can be re-armed without affecting the timing on normal mode channels by pressing function **RUN/STOP**.

### Setting Pulse Timing Parameters

Pulses are defined by a delay, from their sync or start pulse to the active edge, and a width.

- Wid: Sets the width of the active portion of the pulse.
- Dly: Sets the delay from the sync source to the start of the pulse.

*NOTE: If  $Wid + Dly + 75ns$  (hardware reset time)  $> T_o$  Period, the correct pulse width will be generated but at a slower rate.*

### Setting Pulse Output Parameters

There are three basic types of outputs available on the 9520: (a) TTL/CMOS compatible outputs; (b) adjustable amplitude outputs; (c) optical outputs.

Out: Selects between TTL/CMOS mode and Adjustable mode when both are available on a single output.

Pol: Sets the voltage polarity of the pulse, active high or active low. Note: All outputs are positive - negative voltages are not supported.

Ampl: In adjustable mode, it sets the unloaded output voltage. The actual output voltage will depend on the load impedance. For example: If the load is 50 ohms, the output will be 50% of the stated voltage.

### Using the Output Multiplexer

Each output channel includes a multiplexer which allows routing any or all of the timer outputs to the physical output. This allows double pulses and other complex pulse trains to be generated. Only timing parameters are multiplexed together, not amplitudes.

-HGFEDCBA-  
**Mux:** -00000101-

The multiplexer is represented by a “n” bit binary number as shown above. “n” is the number of channels. Each bit represents a channel timer, which is enabled by setting the bit to one. In the above example, timers A and C are combined on the current output.

### Setting System Internal Rate Parameters

The internal  $T_o$  period controls the fundamental output frequency of the system. Each channel may operate at submultiples of the fundamental frequency using their duty cycle mode.

Source: Sets the reference source for the internal  $T_o$  Period.  
 Per: Sets the internal  $T_o$  Period.

To set the system Internal Rate - press the yellow **FUNC** key, then press the **RATE** key, and then use the dial or number pad to specify the  $T_o$  Period.



# 9520 Menus

## 9520 Menu Structure

### System MODE Menus

To Mode	To Mode	To Mode	To Mode
MODE: Continuous	MODE: Single Shot	MODE: Burst	MODE: Duty Cycle
			# / On
		# / Burst	# / Off

### CHANNEL Menus

Timing Parameters	Pulse Configuration	Pulse Configuration
Channel Enable	Channel Enable	Channel Enable
Sync Source	Polarity	Polarity
Pulsewidth	Output Type: TTL, Optical	Output Type: Adjustable
Delay		Output Level

Channel Mode	Channel Mode	Channel Mode	Channel Mode	Wait Menu
Channel Enable	Channel Enable	Channel Enable	Channel Enable	
Mode: Normal	Mode: Single Shot	Mode: Burst	Mode: Duty Cycle	
			# / On	
			# / Off	Wait Count

### Advanced CHANNEL Menus

Multiplexer Menu	Channel Gated Operation	Channel Gated Operation
Channel Enable	Channel Enable	Channel Enable
Mux	Gate Enable	Gate Enable: <method>
Help Line		Gate

RATE Menu

Rate Source	Reference Out
Source	Ref Out
To Period	

TRIG Menu

Trig Mode	Trig Mode
Trig Disabled	
	Threshold
	Edge

GATE Menu

Gate Mode	Gate Mode
Gate Disabled	Gate
	Level

SYSTEM Menus

Comm. Parameters	Comm. Parameters	Comm. Parameters	Keypad Parameters	Misc. Parameters
Interface: RS232	Interface: USB	Interface: GPIB/Ethernet	Key Repeat Rate	Auto Start
Baud Rate		Address	Key Volume	Decimal Mark
Echo Enable			Knob Volume	LCD Brightness

STORE Menu

Store Menu
Configuration #
Name
Help Line

COUNTER Menu

Counter Menu
Enabled/Disabled
T Counts
Help Line

RECALL Menu

Recall Menu
Configuration #
Name
Help Line

INFORMATION Menu

Information Menu
Model Number
Serial Number
Firmware Version
FPGA Version

## System Mode Menu

To Mode	To Mode	To Mode	To Mode
MODE: Continuous	MODE: Single Shot	MODE: Burst	MODE: Duty Cycle
		# / Burst	

### Setting System Mode of Operation

The MODE menu sets the  $T_o$  system timer mode. The menu will show the extra set parameters (Burst, On & Off) only when they are appropriate.

- Mode: Selects the  $T_o$  mode: Continuous, Single Shot, Burst or Duty Cycle mode.
- Burst: Sets the number of pulses to be generated when in Burst mode.
- On: Sets the number of pulses to be generated during each on cycle.
- Off: Sets the number of pulses to skip each during off cycle when in the Duty Cycle mode.

*\*NOTE: Any mode may be started by either the RUN/STOP key in the internal trigger mode or armed by the RUN/STOP key and started by an external trigger in the external trigger mode. In the single shot and burst modes, (internally triggered) the unit disarms itself at the end of the pulse train. Pressing the RUN/STOP key after the unit has been disarmed will generate a new pulse train.*

### Channel Menus

Timing Parameters	Pulse Configuration	Pulse Configuration
Channel Enable	Channel Enable	Channel Enable
Sync Source	Polarity	Polarity
Pulsewidth	Output Type: TTL, Optical	Output Type: Adjustable
Delay		Output Level

Channel Mode	Channel Mode	Channel Mode	Channel Mode	Wait Menu
Channel Enable	Channel Enable	Channel Enable	Channel Enable	
Mode: Normal	Mode: Single Shot	Mode: Burst	Mode: Duty Cycle	
			# / On	
		# / Burst	# / Off	Wait Count

### Enabling Channel Output

At the top of each of the channel menu pages is a parameter to enable or disable the channel. Each channel may be individually controlled. When enabled, the channel key will illuminate.

### Setting the Channel Timing Parameters

To define a pulse requires two parameters: the delay to the active edge and the width of the pulse.

Wid:	Sets the channel pulse width.
Dly:	Sets the channel delay until the active edge.

### Setting Pulse Output Configuration

The 9520 supports two types of outputs: a high speed TTL/CMOS compatible output and for applications which require different voltage levels or higher current, an adjustable voltage output. The pulses can also be defined to be active high or active low.

Pol:	Sets the pulse polarity, active high or active low.
Out:	Selects the output mode, TTL/CMOS or Adjustable
Ampl:	Sets the output voltage level when in the Adjustable mode.

### Setting Channel of Operation

Each channel may be set independently to operate in one of four modes: normal, single shot, burst, or duty cycle (within the **CHANNEL** menus):

Mode:	Selects the mode for the current channel. Additional parameters are provided for the burst mode and the duty cycle mode.
Brst:	Sets the number of pulses in the burst mode to generate before inhibiting output.
On:	Sets the number of pulses to generate before inhibiting output in Duty Cycle Mode.
Off:	Sets the number of pulses to inhibit before repeating the On Cycle in Duty Cycle Mode.

### Delaying the Start of Channel Output

Within any channel mode, the output of the channel can be delayed using the wait parameter (within the **CHANNEL** menu):

Wait:	Sets the number of $T_o$ pulses to wait until enabling the channel output.
-------	--

### Configuring the Channel Multiplexer

To define which channels are fed into the channel multiplexer, the corresponding bit for the desired channel to add should be set to 1.

All desired omitted channels should have the corresponding bit set to 0.

Mux: Enable/disable bitfield.

## Advanced Channel Menus

Advanced Timing Parameters	Advanced Timing Parameters	Channel Gated Operation	Channel Gated Operation
Channel Enable	Channel Enable	Channel Enable	Channel Enable
SyncSource	Sync Source	Gate Enable: Disabled	Gate Enable: <method>
			Gate Polarity

### Setting the Sync Source

Although each channel receives its start pulse from the internal  $T_0$  pulse, the start pulse can be assigned such that the delay entered is relative to the  $T_0$  pulse or any other channel pulse. This allows dependent events to link. The unit will not allow a circular chain of sync sources that would result in a channel triggering itself. The delay entered is relative to the selected sync source.

Sync Source: Selects the channel sync source.

### Setting Channel Gate Control

When the global gate is set (Chan Menu), the channel can then use the gate input with independent behavior from other channels.

Gate: Enables the **GATE** input for the channel by setting the method of output control used with the gating function.  
 Logic: Sets the logic level used with the gating function, either active high or active low.

“Pulse Inhibit” method - the gate prevents the channel from being triggered by the channel’s trigger source pulse. If a pulse has already started when the gate disables the channel, the pulse will continue normal output but will not restart on the next trigger pulse.

“Output Inhibit” method - the gate leaves the base triggering alone and enables/disables the output directly.

### Gate Menu

Gate Mode	Gate Mode
Gate: Disabled	Gate: <method>
	Level
	Logic

### Enabling System Gate

Enables the use of the GATE input as a trigger inhibit or output control for all channels simultaneously, or on a per channel basis.

Mode: Selects between disabling the **GATE** inputs and method of output control  
Level: Sets the gating threshold.  
Logic: Sets the active logic level.

### Trig Menu

Trig Mode	Trig Mode
Trig: Disabled	Trig: <trig mode>
	Level
	Edge

### Enabling System Trigger

Enable the use of the **TRIG** input by the system timer as a trigger source.

Mode: Selects between disabling/enabling the trigger mode(s).  
Level: Sets the trigger threshold.  
Edge: Selects between rising edge / falling edge as the trigger source when a trigger mode is enabled.

### Rate Menu

Rate Source	Reference Out
Source	Ref Out
To Period	

### Setting the Clock Source and Internal Rate

Source: Selects the internal or external clock source from which the unit will operate.  
Per: Sets the  $T_o$  period which determines the fundamental output frequency of the unit.

### Setting the Output Reference

Ref Out: Selects the frequency of the output reference for synchronizing with external system components.

## System Menu

Comm. Parameters	Comm. Parameters	Comm. Parameters	Keypad Parameters	Misc. Parameters
Interface: RS232	Interface: USB	Interface: GPIB/Ethernet	Key Repeat Rate	Auto Start
Baud Rate		Address	Key Volume	Decimal Mark
Echo Enable			Knob Volume	LCD Brightness

### Setting System Communication Parameters

The 9520 comes with a standard RS232 serial port and USB port. The unit will not respond to computer commands unless these ports are properly configured.

Interface:	RS232, USB, GPIB (optional), Ethernet (optional)
Baud Rate:	Selects the baud rate for the selected interface.
Echo:	Selects whether to echo characters back to the host computer or not.
Address:	Sets the GPIB address.

### Setting Keypad Parameters

The rate at which a key will repeat itself when held down may be set. This can be used to provide a controlled rate at which a parameter is incremented. In addition, the volume of the beep can be controlled for both the keypad and the adjustable knob.

Key Rate:	Sets the rate at which the keys will repeat when held down.
Key Vol:	Sets the beep volume for the keypad.
Knob Vol:	Sets the beep volume for the Rotary Knob.

### Setting the Auto Start Mode

The unit may be configured to automatically start generating pulses after power up.

### Setting the Display Decimal Mark

Mark:	Selects the format of the decimal mark, "." or ",".
-------	---

### Setting the Display Brightness

LCD	Adjusts display brightness.
-----	-----------------------------

## Store Menu

<b>Store Menu</b>
Configuration #
Name
Help Line

### Storing a Configuration

Use the following procedure to store a complete system configuration:

- Set all parameters to the desired value.
- Select a configuration number.

*\*NOTE: You cannot store to the zero location, as that contains the factory default values.*

- Label the configuration as desired.
- From the Store menu, press the store button sequence (FUNC + store).

## Recall Menu

<b>Recall Menu</b>
Configuration #
Name
Help Line

### Recalling System Configurations

Use the following procedure to recall a stored or default system configuration:

- Enter the Recall Menu (FUNC + recall).
- Select a configuration number.
- From the Recall Menu, press the recall key sequence (function + recall).

*\*Note: Configuration 0 is the factory default setting.*

## Counter Menu

<b>Counter Menu</b>
Enable/Disable
T Counts
Help Line

### Using the Counter Function

The Counter function counts the number of  $T_o$  pulses output by the system clock. When the unit is operated in system single shot mode, the  $T_o$  count reflects the number of incoming trigger pulses.

## Information Menu

<b>Information Menu</b>
Model Number
Serial Number
Firmware Version
FPGA Version

The Information Menu provides all of the pertinent version numbers and serial numbers for the unit. This information should be readily available when contacting customer service for troubleshooting help.



---

# Operating the 9520

## Quick Start - Normal Internal Rate Generator Operation

The 9520 has a powerful set of functions providing a number of modes of operation for the internal or “System” rate generator ( $T_o$ ). Most of these functions can be ignored if a simple continuous stream of pulses is required. Starting from the default settings, which can be restored by recalling configuration 0, the following parameters need to be set:

Pulse Width, Delay	Enter the Channel menus by pressing the letter key. Enter the required pulse width and delay. Repeat for each output channel.
$T_o$ Period	Enter the Rate menu by pressing the <b>FUNC</b> key and then the <b>RATE</b> key. Set the desired pulse period. Note that in general, the pulse delay plus the pulse width, plus a 75ns hardware reset constant, for any channel must be less than the $T_o$ period.
Start	Press the <b>RUN/STOP</b> key to start generating pulses.
Stop	Press the <b>RUN/STOP</b> key a second time to stop generating pulses.

## Quick Start - Normal External Trigger Operation

To generate a single pulse for every external trigger event, based on the default configuration 0, the following parameters need to be set:

System Mode	Enter the <b>MODE</b> menu by pressing the <b>FUNC</b> key and then the <b>MODE</b> key. Select Single Shot mode.
Trig	Enter the <b>TRIG</b> menu by pressing the <b>TRIG</b> key. Select Triggered.
Level	Press the <b>NEXT</b> key until the Level parameter is highlighted. Set the trigger threshold voltage to approximately 50% of the trigger signal amplitude.

Edge	Press the <b>NEXT</b> key until the Edge parameter is highlighted. Set the instrument to trigger off the rising edge or falling edge as desired.
Pulse Width, Delay	Enter the Channel menus by pressing the letter key. Enter the required pulse width and delay. Repeat for each output channel.
Start	Press the <b>RUN/STOP</b> key to start/arm the instrument. The 9520 will now generate a pulse for every valid trigger.
Stop	Press the <b>RUN/STOP</b> key a second time to stop/disarm the instrument (i.e. to stop generating pulses).

## System Timer Overview

For internal operation, the 9520 contains a timer and mode generator which generates an internal  $T_0$  clock that is used to trigger all the channel timers. System modes are controlled via the **MODE** menu.

### To Use Continuous Mode

The **RUN/STOP** button starts and stops a continuous pulse stream at the rate specified by the Rate menu. This corresponds to the normal output mode for most pulse generators.

To generate a continuous stream of pulses:

- within the system **Mode** menu

Mode                                      Select Continuous for the system mode.

- within the **Rate** menu

Source                                      Select the system oscillator or the external clock in frequency.

Period                                      Set the desired period.

Pressing the **RUN/STOP** key will now generate a stream of  $T_0$  pulses, at a rate specified by the period parameter.

### To Use Single Shot Mode

To generate a single pulse with every press of the **RUN/STOP** key:

- within the system **Mode** menu

Mode                                      Select Single Shot for the system mode.

Pressing the **RUN/STOP** key will now generate a single pulse.

### To Use System Burst Mode Function

The **RUN/STOP** button generates a stream of “n”  $T_o$  pulses, where “n” is specified by the Burst parameter. The rate is specified in the Rate menu. Pressing the **RUN/STOP** button while the burst is in process will stop the output. After the burst has been completed, pressing the **RUN/STOP** button will generate another burst. To generate a burst of pulses:

- within the system **Mode** menu

Mode: Select the Burst mode.

Burst: Set the number of pulses to produce in the burst.

### To Use System Duty Cycle Function

The **RUN/STOP** button starts a continuous pulse stream which oscillates on for the “n” pulses and off for “m” pulses, where “n” and “m” are specified by the On and Off parameters, respectively. The rate is specified in the Rate Menu. To generate a stream of pulses which oscillates on for ‘n’ pulses and off for ‘m’ pulses:

- within the system **Mode** menu

Mode: Select the Duty Cycle mode.

On: Set the number of pulses to produce during the on cycle.

Off: Set the number of pulses to skip during the off cycle.

-within the **Rate** menu

Source Select the system oscillator or the external clock in frequency.

Period: Set desired Period.

### Channel Timer Overview

The output of each channel is controlled by two timers to generate the delay timing and the pulse width. All channels are simultaneously triggered, depending on the system mode, by either the internal  $T_o$  pulse, the external trigger, or a trigger provided by a cpu. A given channel may or may not generate a pulse depending on its own channel mode as described below.

When one channel is generating a continuous stream of pulses, a user can trigger a single shot or burst of pulses on another channel without interrupting the continuous stream by pressing “**FUNC**” and “**RUN/STOP**”.

### To Use Channel Normal Mode Function

The Normal mode generates a continuous stream of pulses at a rate determined by the system timer:

- within the **Channel** menus

Enable: Select Enable to enable channel output.

Dly: Set the desired delay.

Wid: Set the desired pulse width.  
Mode: Select the Normal mode.

Pressing the **RUN/STOP** key will now generate a continuous stream of pulses.

### To Use Channel Single Shot Function

The Single Shot mode generates a single pulse every time the **RUN/STOP** key is pressed. If the unit is in the active state, ( i.e. channels which are set to the Normal mode are producing pulses), pressing the “**FUNC**” key and **RUN/STOP** key will reset the Single Shot counters and generate one pulse in sync with the other channels running in the Normal mode. To use the Single Shot mode:

- within the **Channel** menus

Enable Select Enable to enable channel output.  
Delay Set the desired delay.  
Width Set the desired pulse width.  
Mode Select the Single shot mode.

### To Use Channel Burst Mode Function

The Burst mode generates a burst of pulses every time the **RUN/STOP** key is pressed. If the unit is in the active state, (i.e. channels which are set to the Normal mode are producing pulses), pressing the “**FUNC- RUN/STOP**” key sequence will reset the Burst counters and generate a new set of pulses in sync with the other channels running in the Normal mode. “**FUNC - RUN/STOP**” will not affect  $T_o$  pulse status. To use the Burst mode:

-within the **Channel** menus

Enable Select Enable to enable channel output.  
Delay Set the desired delay.  
Width Set the desired pulse width.  
Mode Select the Burst mode.  
#/Burst Set the number of pulses to produce in the burst.

### To Use the Channel Duty Cycle Function

To generate a stream of pulses which oscillates on for ‘n’pulses and off for ‘m’ pulses:

- within the **Channel** menus

Enable Select Enable to enable channel output.  
Delay Set the desired delay.  
Width Set the desired pulse width.  
Mode Select the Duty Cycle mode.  
On Cycle Set the number of pulses to produce during the on cycle.  
Off Cycle Set the number of pulses to skip during the off cycle.

*Note: Older Quantum Composers pulse generators had a divide-by-n function. The duty cycle mode is a more general case. To reproduce the divide-by-n function, set the on cycle to 1 and set the off cycle to (n-1), where “n” is the divide-by-n factor.*

### **To Use the Channel Gating Function**

Each channel may use the external input to gate or control its output. The gate controls the triggering of the channel. To use the gate, set the following parameters

- within the **Channel** menu

Mode: In Gate Menu must be set to Channel Menu.  
Channel Gate: Select Pulse Inh or Output Inh.  
Logic: Select active high or active low.

In “Pulse” inhibit method, the gate prevents the channel from being triggered by the channel’s trigger source pulse. If a pulse has already started when the gate disables the channel, the pulse will continue normal output but will not restart on the next pulse. In “Output” inhibit method, the gate leaves the base triggering alone and enables/disables the output directly. Output pulses will immediately cease when the gate signal is removed

### **External Input Overview**

The external inputs may be used to trigger the unit or to gate the system or channel timers. When using a trigger input, the external input acts as a system start pulse. Depending on the system mode, the result of a trigger input can be either a single pulse, a burst of pulses or the start of a stream of pulses.

### **To Generate a Pulse on Every Trigger Input**

To generate a pulse on every external trigger received, set the following parameters:

- within the **Mode** menu

Mode: Select the Single Shot mode.

- within the **Trigger** menu

Mode: Select Triggered mode.  
Level: Set the trigger threshold level.  
Edge: Select which edge, rising or falling, to trigger on.

Pressing the **RUN/STOP** key will arm the unit. Once the unit is armed, it will generate a  $T_0$  pulse for every external trigger received. Pressing the **RUN/STOP** key will disarm the unit. This mode corresponds to the normal external trigger mode found on most other pulse generators.

### To Generate a Burst of Pulses on Every Trigger Input

To generate a burst of pulses for every external trigger received set the following parameters:

-within the **Mode** menu

- Mode: Select the Burst mode.
- Burst: Set the number of pulses to generate in each burst.-  
within the Rate menu
- $T_o$ : Set the period between pulses.

- within the **Rate** menu

- Source: Select the system oscillator or an external clock in frequency.
- $T_o$ : Set the period between pulses.

- within the **Trigger** menu

- Mode: Select Triggered mode.
- Level: Set the trigger threshold level.
- Edge: Select which edge, rising or falling, to trigger on.

Pressing the **RUN/STOP** key will arm the unit. Once the unit is armed, it will generate a set of pulses for every external trigger received. The unit is reset at the end of a burst and will generate another set of pulses upon receiving a new trigger. Triggers that occur in the middle of a burst are ignored. Pressing the **RUN/STOP** key will disarm the unit.

### To Start a Continuous Stream of Pulses Using the External Trigger

The external trigger may be used to start the unit generating pulses:

- within the **Mode** menu

- Mode: Select the Continuous mode.

- within the **Rate** menu

- Source: Select the system oscillator or an external clock in frequency.
- $T_o$ : Set the period between pulses.

- within the **Trigger** menu

- Mode: Select the Trigger mode.
- Level: Set the trigger threshold level.
- Edge: Select which edge, rising or falling, to trigger on.

Pressing the **RUN/STOP** key will arm the unit. Once the unit is armed, it will begin generating pulses after an external trigger is received. Triggers that occur after the pulses start are ignored. Pressing the **RUN/STOP** key a second time will disarm the unit.

### To use the External Gate to Control the System

The external gate may be used to control the output of the unit. To gate the system timer:

- within the **Mode** menu

Mode: Select the desired mode.

- within the **Rate** menu

Source: Select the system oscillator or an external clock in frequency..

Per: Set the period between pulses.

- within the **Gate** menu

Mode: Select Pulse Inh or Output Inh.

Level: Set the gate threshold level.

Logic: Select active high or active low.

Pressing the **RUN/STOP** key will arm the unit. Once the unit is armed, it will begin generating pulses whenever the external gate input is in the active state. Pressing the **RUN/STOP** key a second time will disarm the unit.



---

# Programming the 9520

## Personal Computer to Pulse Generator Communication

The 9520 comes standard with an RS232 serial and USB interface. An Ethernet and GPIB interface is available as an option. All menu settings can be set and retrieved over the computer interface using a simple command language. The command set is structured to be consistent with the Standard Commands for Programmable Instruments. Although due to the high number of special features found in the 9520, many of the commands are not included in the specification. The syntax is the same for all interfaces.

## RS232 Interface Overview

The serial port is located on the back of the 9520 and uses a 9-pin D-type connector with the following pinout (as viewed from the back of the unit):

1	No Connection
2	Tx - Transmit (to computer)
3	Rx - Receive (from computer)
4	DTR - Connected to pin 6
5	Ground
6	DSR - Connected to pin 4
7	RTS - Connected to pin 8
8	CTS - Connected to pin 7
9	No Connection

The serial port parameters should be set as follows:

Baud Rate	4800, 9600 19200, 38400, 57600, 115200*
Data Bits	8
Parity	None
Stop Bits	1

\*The default baud rate for the RS232 is 115200.

## USB Interface Overview

The USB interface is standard on the 9520. Before this type of communication can be used, the appropriate drivers must be installed on the personal computer (pc). These drivers are included on the CD that was shipped with your unit. Please contact Quantum Composers or visit [www.quantumcomposers.com](http://www.quantumcomposers.com) for updated installation files and instructions.

USB communication is achieved by using a mapped (virtual) COM port on the PC. The driver installation executable will obtain an unused COM port number,

install the USB drivers, and make that COM port number available for typical RS232 communication to the pulse generator. Hyperterminal or other common software may be used.

When communicating through the mapped COM port over USB, the baud rate for the communication port used by the USB chip must match the baud rate for the COM port on the PC. Access to the USB port baud rate is done using the SCPI command “:SYSTem:COMMunicate:SERial:USB n” command, where ‘n’ is the desired communication speed. This parameter can be accessed via any communication method. The default baud rate for USB is 38400.

USB communication notes:

- The correct drivers must be installed on the personal computer before communication can be accomplished via USB.
- The BAUD rates on the PC and on the pulse generator must match for successful communication.
- The USB port’s BAUD rate on the pulse generator can be set using the SCPI command “:SYSTem:COMMunicate:SERial:USB n” where ‘n’ can be:
  - o 4800
  - o 9600
  - o 19200
  - o 38400 (default)
  - o 57600
  - o 115200
- USB 1.0 specification is used. The USB cable can be removed without “unplugging” the device in the operating system environment.
- Echo functionality is not available on the USB port.

## **GPIB Interface Overview**

Also known as IEEE-488, a GPIB computer interface is optional on the 9520. Before using this interface, the address must be set using the GPIB address menu item. The command set is the same for the RS-232, GPIB, USB and Ethernet. Different interfaces may be used at the same time. Responses will be made to the most recently used interface. Refer to Appendix F included at the end of this manual for more information about the GPIB Interface and Operation.

## **Ethernet Interface Overview**

An Ethernet interface is optional on the 9520. Refer to Appendix G included at the end of this manual for more information about the Ethernet Interface and Operation.

## Programming Command Types and Format

The 9520 Pulse Generator uses two types of programming commands: IEEE 488.2 Common Commands and Standard Commands for Programmable Instruments (SCPI). The format is the same for all interfaces. Hyperterminal (in Windows) or any other generic terminal program may be used to interactively test the commands using the RS232 interface. The format of each type is described in the following paragraphs.

### Line Termination

The pulse generator uses text-style line terminations. When a command is sent to the unit, the firmware is programmed to read characters from a communication port until it reads the line termination sequence.

The command string is parsed and executed after reading these characters. These characters are the “carriage return” and “linefeed”. They are ASCII character set values 13 and 10 respectively (hex 0x0D and 0x0A). All command strings need to have these characters appended.

When the pulse generator responds to a command, whether it is a query or a parameter change, it also appends its return strings with these characters. Coded applications could use this behavior to know when to stop reading from the unit. However, if the “echo” parameter is enabled, there will be two sets of line terminators, one following the echoed command string, and one following the pulse generator’s response.

*Note: The pulse generator will echo commands on the DB9 serial port only.*

The pulse generator responds to every communication string. If the communication string is a query, the unit responds with the queried response (or error code) followed by the line terminators. If the communication string is a parameter change, the response is “ok” (or error code) followed by the line terminators. For this reason, it is not recommended that multiple commands be stacked together into single strings as is common with some other types of instruments. It is recommended that the coded application send a single command in a string and follow immediately by reading the response from the unit. Repeat this sequence for multiple commands.

### IEEE 488.2 Common Command Format

The IEEE 488.2 Common Commands control and manage generic system functions such as reset, configuration storage and identification. Common commands always begin with the asterisk (\*) character and may include parameters. The parameters are separated from the command mnemonic by a space character. For Example:

```
*RST      <cr> <lf>
*RCL 1    <cr> <lf>
*IDN?     <cr> <lf>
```

## SCPI Command Keywords

The commands are shown as a mixture of upper and lower case letters. The upper case letters indicate the abbreviated spelling for the command. You may send either the abbreviated version or the entire keyword. Upper and/or lower case characters are acceptable.

For example, if the command keyword is given as POLarity, then POL and POLARITY are both acceptable forms; truncated forms such as POLAR will generate an error; polarity, pol, and PolAriTy are all acceptable as the pulse generator is not case sensitive.

## SCPI Command Format

SCPI commands control and set instrument specific functions such as setting the pulse width, delay and period. SCPI commands have a hierarchical structure composed of functional elements that include a header or keywords separated with a colon, data parameters, and terminators. For example:

### SCPI Format

```
:PULSE1:STATE ON           <cr> <lf>
:PULSe1:WIDTh 0.000120     <cr> <lf>
:PULSe:POL NORMal         <cr> <lf>
```

Any parameter may be queried by sending the command with a question mark appended. For example:

### - QUERY FORMAT

```
:PULSE1:STATE?           <cr><lf>
    Will return:         <cr><lf>

:PULSE1:WIDT?           <cr><lf>
    Will return:         0.000120000 <cr><lf>

:PULSE1:POL?           <cr><lf>
    Will return:         NORM <cr><lf>
```

## SCPI Keyword Separator

A colon (:) must always separate one keyword from the next lower-level keyword. A space must be used to separate the keyword header from the first parameter. If more than one parameter is used, you must separate subsequent parameters with a comma.

## SCPI Optional Keywords

Optional keywords and/or parameters appear in square brackets ( [ ] ) in the command syntax. Note that the brackets are not part of the command and should not be sent to the pulse generator. When sending a second level keyword without the optional keyword, the pulse generator assumes that you intend to use the optional keyword and responds as if it had been sent.

## SCPI Specific and Implied Channel

Some commands, such as PULSe, allow specifying a channel with an optional numeric keyword suffix. The suffix will be shown in square brackets [ 1 / 2 ]. The brackets are not part of command and are not to be sent to the pulse generator. The numeric parameters correspond to the following channels: 0 = To, 1 = ChA, 2 = ChB, etc. Only one channel may be specified at a time.

If you do not specify the channel number, the implied channel is specified by the :INSTrument:SELEct command or the last referenced channel.

After power-up or reset (\*RST) The instrument default is channel #1.

## SCPI Parameter Types

The following parameter types are used:

<numeric value>	Accepts all commonly used decimal representation of numbers including optional signs, decimal points, and scientific notation: 123, 123e2, -123, -1.23e2, .123, 1.23e-2, 1.2300E-01.
<boolean value>	Represents a single binary condition that is either true or false. True is represented by a 1 or ON; false is represented by a 0 or OFF. Queries return 1 or 0.
<identifier>	Selects from a finite number of predefined strings.

## Error Codes

The unit responds to all commands with either:  
ok <cr><lf> or ?n <cr><lf>

Where "n" is one of the following error codes:

- 1 Incorrect prefix, i.e. no colon or \* to start command.
- 2 Missing command keyword.
- 3 Invalid command keyword.
- 4 Missing parameter.
- 5 Invalid parameter.
- 6 Query only, command needs a question mark.
- 7 Invalid query, command does not have a query form.
- 8 Command unavailable in current system state.

## Programming Examples

Example 1) 20 ms pulse width, 2.3 ms delay, 10 Hz, internal trigger, continuous operation.

:PULSE1:STATE ON <cr> <lf>	enables channel A
:PULSE1:POL NORM <cr> <lf>	sets polarity to active high
:PULSE:WIDT 0.020 <cr> <lf>	sets pulse width to 20 ms
:PULSE1:DELAY 0.0023 <cr> <lf>	sets delay to 2.3 ms
:PULSE0:MODE NORM <cr> <lf>	sets system mode to continuous
:PULSE0:PER 0.1 <cr> <lf>	sets period to 100 ms (10 Hz)
:PULSE0:TRIG:MODE DIS <cr> <lf>	disables the external trigger

To start the pulses use either of the following commands:

:PULSE0:STATE ON <cr> <lf>	starts the pulses
:INST:STATE ON <cr> <lf>	alternate form to start pulses

Example 2) 25 $\mu$ s pulse width, 0 delay, external trigger, one pulse for every trigger.

:PULSE1:STATE ON <cr> <lf>	enables channel A
:PULSE1:POL NORM <cr> <lf>	sets polarity to active high
:PULSE:WIDT 0.000025 <cr> <lf>	sets pulse width to 25 $\mu$ s
:PULSE1:DELAY 0 <cr> <lf>	sets delay to 0
:PULSE0:MODE SING <cr> <lf>	sets system mode to single shot
:PULSE:TRIG:MODE TRIG <cr> <lf>	sets system to external trigger
:PULS:TRIG:LEV 2.5 <cr> <lf>	sets trigger level to 2.5 v
:PULS:TRIG:EDGE RIS <cr> <lf>	set to trigger on rising edge

To arm the instrument in external gate mode, use either of the following commands:

:PULSE0:STATE ON <cr> <lf>	arms the instrument
:INST:STATE ON <cr> <lf>	alternate form if T <sub>0</sub> is currently selected

A software generated external trigger can be generated by using the following command:

*TRG <cr> <lf>	generates a software external trigger
----------------	---------------------------------------

## 9520 SCPI Command Summary

Keyword	Parameter	Std/ New	Comments
:INSTrument		Std	Subsystem. Supports treating each channel as a logical instrument.
:CATalog?		Std	Query only. Returns a comma-separated list of the names of all channels. A two channel instrument would return: T <sub>0</sub> , CHA, CHB.
:FULL?		Std	Query only. Returns a comma-separated list of the names of all channels and their associated number. A two channel instrument would return: T <sub>0</sub> , 0, CHA, 1, CHB, 2.
:COMMands?		New	Query only. Returns an indented list of all SCPI commands.
:NSElect	<numeric value>	Std	Selects a channel using the channel's numeric value. All channel specific commands will refer to the selected channel.
:SElect	<identifier>	Std	Selects a channel using the channel's identifier string. All subsequent channel specific commands will refer to the selected channel.
:STATe	<boolean value>	Std	Enables/Disables the selected channel output. If T <sub>0</sub> is selected all output is affected. Enabling T <sub>0</sub> is the same as pressing the RUN button.

## 9520 SCPI Command Summary

Keyword	Parameter	Std/ New	Comments
:PULSe[0]		Std	Subsystem. Contains commands to control the output pulse generation. Commands without suffix refer to the currently selected logical instrument. See INSTrument subsystem.
:COUNter		New	Subsystem. Contains commands to define the Counter function.
:STATe	<boolean value>	New	Enables/Disables the counter function.
:CLear	TCNTS/GCNTS	New	Clears the designated counter. Standard units only have the Trigger counter.
:COUNter?	TCNTS/GCNTS	New	Queries the number of counts for the specified input. Standard units only have the Trigger counter.

## 9520 SCPI Command Summary

Keyword	Parameter	Std/ New	Comments
:STATe	<boolean value>	Std	Enables / Disables the output for all channels. Command is the same as pressing the RUN/STOP button.
:PERiod	<numeric value>	Std	Sets the To period.
:MODE	NORMAL / SINGLE / BURSt / DCYClE	New	Sets the To mode.
:BCOunter	<numeric value>	New	Burst Counter. Number of pulses to generate in the Burst mode.
:PCOunter	<numeric value>	New	Pulse Counter. Number of pulses to generate during on cycle of the Duty Cycle mode.
:OCOunter	<numeric value>	New	Off Counter. Number of pulses to inhibit output during the off cycle of the Duty Cycle mode.
:ICLock	SYS/ EXT10/ EXT20/ EXT25/ EXT40/ EXT50/ EXT80/ EXT100	New	Sets Source for the internal rate generator. System Clock or External Source ranging from 10MHz to 100MHz.
:OCLock	To / 10 / 11 / 12 / 14 / 16 / 20 / 25 / 33 / 50 / 100	New	Sets external clock output. To Pulse or 50% duty cycle TTL output from 10MHz to 100MHz.
:GATe		New	Subsystem. Contains the commands to define the Gate function.
:MODE	DIS/PULS/OUTP/ CHAN	New	Sets Global Gate Mode. Disable, pulse inhibit, output inhibit, channel.
:LOGic	LOW / HIGH	New	Sets Channel Gate logic level. Active low or active high.
:LEVel	<numeric value>	New	Sets the gate threshold. Value is in volts with a range of .20 to 15 Volts.
:TRIGger		New	Subsystem. Contains the commands to define the Trigger function.
:MODE	DIS / TRIG	New	Sets Trigger Mode. Disable or TRIG (enable).
:EDGE	RISing / FALLing	New	Selects which edge (rising or falling) to use as the trigger signal.
:LEVel	<numeric value>	New	Sets the Trigger Threshold. Value is in volts, with a a range of .20 to 15 Volts.

## 9520 SCPI Command Summary

Keyword	Parameter	Std/ New	Comments
:PULSe [1 / 2 / n]		Std	Subsystem. Contains commands to control the output pulse generation. Valid suffix range depends on the number of channels (ChA = 1, ChB = 2, etc). Command without suffix refers to the currently selected logical instrument. See INSTRUMENT subsystem.
:STATe	<boolean value>	Std	Enables/Disables the output pulse for selected channel.
:WIDTh	<numeric value>	Std	Sets the width or duration of the output pulse.
:DELay	<numeric value>	Std	Sets the time from the start of the To period to the first edge of the pulse.
:SYNC	TO, CHA, CHB, CHC, CHD, etc.	New	Selects the Sync source.
:MUX	<numeric value>	New	Selects which timers are enabled as output for the current channel.
:POLarity	NORMAL / COMPLEMENT / INVERTed	Std	Sets the polarity of the pulse. For NORMAL operation the second nominal state is more positive than the first. COMPLEMENT and INVERTed are aliases. For both, the second state is more negative than the first.
:OUTPut		New	Subsystem. Contains command to control output mode.
:MODE	TTL / ADJUSTable /	New	Selects output AMPLitude mode: TTL/CMOS, ADJUSTable.
:AMP	<numeric value>	New	Sets adjustable output level.
:CMODE	NORMAL / SINGLE / BURSt / DCYCLE	New	Channel Mode. Sets the channel pulse series output mode.
:BCOUNTER	<numeric value>	New	Burst Counter. Sets the number of pulses to generate when channel is in the BURST mode.
:PCOUNTER	<numeric value>	New	Pulse Counter. Sets the number of pulses to generate during the on cycle of the Duty Cycle Mode.
:OCOUNTER	<numeric value>	New	Off Counter. Number of pulses to inhibit output during the off cycle of the Duty Cycle mode.
:WCOUNTER	<numeric value>	New	Sets the number of To pulses to delay until enabling output.
:CGATE	DIS / PULS / OUTP	New	Sets Channel Gate Mode. Disable, pulse inhibit, output inhibit. (Global Gate Mode must be set to CHAN for this command to be available).
:CLOGic	LOW / HIGH	New	Sets Channel Gate Logic level. Active low or active high. (Global Gate Mode must be set to CHAN for this command to be available).

## 9520 SCPI Command Summary

Keyword	Parameter	Std/ New	Comments
:SYSTem		Std	
:STATe?		New	Query only. Returns the state of the machine: returns "1" if the machine is armed and/or generating pulses or "0" if the machine has been disarmed.
:BEEPer		Std	Subsystem. Controls the audible beeper.
:STATe	<boolean value>	Std	Enables/disables the beeper.
:VOLume	<numeric value>	Std	Sets the volume of the beeper. Range is 0 to 100, where 0 is off and 100 is maximum volume.
:COMMunicate		Std	Subsystem. Controls the RS232 and GPIB interfaces.
:GPIB		Std	Subsystem. Controls the physical configuration of the GPIB port.
:ADDRes	<numeric value>	Std	Sets the GPIB of the instrument.
:SERial		Std	Subsystem. Controls the physical configuration of the RS232 port.
:BAUD	4800 / 9600 / 19200 / 38400 / 57600 / 115200 /	Std	Sets the baud rate for both receiving and transmitting using the DB9 RS232 port.
:USB	4800 / 9600 / 19200 / 38400 / 57600 / 115200 /	New	Sets the baud rate for communication when using mapped comports for USB and Ethernet communication. Must be set to the default value (115200) for raw TCP/IP communication (i.e. Labview).
:ECHO	<boolean value>	New	Enables/Disables transmission of characters received on the DB9 serial port.
:KLOCK	<boolean value>	New	Locks the keypad.
:AUTorun	<boolean value>	New	After power-up, unit will start generating pulses automatically.
:VERSion?		Std	Query only. Returns SCPI version number in the form: YYYY.V ex. 1999.0
:SERN?		New	Query only. Returns the serial number.
:INFORmation?		New	Query only. Returns model, serial number, firmware version, and FPGA version numbers.
:NSID?		New	Query only. Returns firmware and FPGA identification numbers.
:CAPS	<boolean value>	New	Forces unit to recognize commands only sent in capital letters. 1 turns on the feature - 0 disables the feature.

Keyword	Parameter	Std/ New	Comments
:DISPlay		Std	Subsystem. Contains commands to control the display.
:MODE	<boolean value>	New	Enables/Disables automatic display update. When true, front panel display is updated with serial command parameter changes. Setting to false decreases response time.
:UPDate?		New	Query only. Forces update of display. Use when mode is false.
:BRiGhtness	<numeric value>	New	Controls intensity of display. Range is 0 to 4, where 0 is off and 4 is full intensity.
:ENABle	<boolean value>	New	Enables/Disables the display and front panel lights. When Disabled the keylock is enabled to prevent parameter changes from the front panel.

## IEEE 488.2 Common Commands

Mnemonic	Command Name	Parameters	Comments
*IDN?	Identification Query		Queries the Pulse Generator Identification. The ID will be in the following format: model#-#channels-option#-version#
*RCL	Recall Command	<numeric value>	Restores the state of the Pulse Generator from a copy stored in local nonvolatile memory (0 through 12 are valid memory blocks).
*RST	Reset Command		Resets the Pulse Generator to the default state.
*SAV	Save Command	<numeric value>	Stores the current state of the Pulse Generator in local nonvolatile memory (1 through 12 are valid memory blocks).
*TRG	Trigger		Generates a software trigger pulse. Operation is the same as receiving an external trigger pulse.
*LBL	Setup Label	<string value>	Query Form returns the label of the last saved or recalled configuration.  Command Form sets the label string for the next "**sav" command. String must be in double quotes, 14 characters max.
*ARM	Channel Trigger Reset		Resets channel triggers when channels are set to single shot or burst mode. Functions like pressing the function then run/stop button.

---

---

## Appendix

---

# Appendix A - Specifications

## 9520 Specifications

### PULSE GENERATION

RANGE	
DELAY	0 - 1000s
WIDTH	10ns - 1000s
RESOLUTION	250ps
TIMEBASE	50MHz
TIMEBASE ACCURACY	25PPM

### INTERNAL RATE GENERATOR

RATE ( $T_o$ period)	0.0002Hz to 10.000MHz
RESOLUTION	10ns
ACCURACY	Same as timebase
JITTER	50ps
BURST MODE	1 to 9,999,999

### TTL/ADJUSTABLE OUTPUTS

NUMBER	2, 4 or 8 Channel Output
LOAD	50 ohm
RISE TIME (10% - 90%)	3ns typ TTL 15ns typ @ 20V (high imp) Adj 25ns typ @ 10V (50 ohms) Adj
SLEW RATE	>0.5V/ns TTL >0.1V/ns Adj
OVERSHOOT	<100mV + 10% of pulse amplitude
LEVELS	TTL 0 to 4 VDC high imp 2.0 to 20.0 VDC high imp 1.0 to 10.0 VDC into 50 ohm
RESOLUTION	10mV
ACCURACY	1ns + .0001 x Setpoint

## **EXTERNAL INPUTS**

### **TRIGGER INPUTS**

NUMBER	0, 2 or 4
RATE	DC to 1/ (200ns + longest delay)
THRESHOLD	0.2 to 15VDC
MAXIMUM INPUT VOLT.	60V Peak
RESOLUTION	10mV
IMPEDANCE	1000 ohm
SLOPE	Rising or Falling
IMPEDANCE	1000 ohm
JITTER	800ps RMS
INSERTION DELAY	<180ns
MINIMUM PULSE WIDTH	2ns

### **GATE INPUTS**

THRESHOLD	0.2 to 15 VDC
MAXIMUM INPUT VOLT.	60V Peak
RESOLUTION	10mV
POLARITY	Active High/Active Low
FUNCTION	Pulse Inhibit or Output Inhibit
CHANNEL BEHAVIOR	Global w/Individual Channel
PULSE INHIBIT DELAY	120ns
OUTPUT INHIBIT DELAY	50ns

## **OPTICAL OUTPUTS**

NUMBER	2, 4 or 8
WAVELENGTH	820nm or 1300nm
MAXIMUM SIGNAL RATE	5MBd
MAXIMUM LINK DIST.	1.5km
CONNECTOR TYPE	ST
ACCURACY	1ns + .0001 x Setpoint

## **OPTICAL INPUTS**

NUMBER	0, 2 or 4
WAVELENGTH	820nm or 1300nm
MAXIMUM SIGNAL RATE	5MBd
MAXIMUM LINK DIST.	1.5km
CONNECTOR TYPE	ST
INSERTION DELAY	<300ns
JITTER	<15ns

## **OUTPUT MODULES**

### Standard

AT20 Dual channel, TTL/CMOS & Adjustable output module

### Optional

L82 Dual channel, 820nm optical output module

L130 Dual channel, 1300nm optical output module

AT35 Dual channel, TTL/35V high voltage output module

TZ50 Dual channel, high current TTL/CMOS (for driving 50 ohm loads) & adjustable output module

TZ35 Dual channel, high current TTL/CMOS (for driving 50 ohm loads) & 35V high voltage output module

DD25 Dual channel, 1.0 A LED constant current (PWM controlled) driver module

## **INPUT MODULES**

### Standard

IA15 Dual channel, 1 trigger / 1 gate input module

### Optional

IL82 Dual channel, 820nm optical input module

IL130 Dual channel, 1300nm optical input module

### System Options

I Incrementing (provides automatic high speed incrementing/ decrementing of delay and/or pulse width for each channel)

DT15 Dual Trigger Logic – provides additional trigger via gate input

COM Extended Communications – Adds Ethernet & GPIB

SRM Single Rackmount

DRM Dual Rackmount

***\*Other custom modules (LED drivers, higher voltages, current sources) available, call with your request.***

# Appendix B - Safety Symbols

## Safety Marking Symbols

Technical specifications including electrical ratings and weight are included within the manual. See the Table of Contents to locate the specifications and other product information. The following classifications are standard across all QC products:

- Indoor use only
- Ordinary Protection: This product is NOT protected against the harmful ingress of moisture.
- Class 1 Equipment (grounded type)
- Main supply voltage fluctuations are not to exceed +/-10% of the nominal supply voltage.
- Pollution Degree 2
- Installation (overvoltage) Category II for transient overvoltages
- Maximum Relative Humidity: <80% RH, non-condensing
- Operating temperature range of 0° C to 40° C
- Storage and transportation temperature of -40° C to 70° C
- Maximum altitude: 3000 m (9843 ft.)
- This equipment is suitable for continuous operation.

This section provides a description of the safety marking symbols that appear on the instrument. These symbols provide information about potentially dangerous situations which can result in death, injury, or damage to the instrument and other components.

**Symbols      Publications; Descriptions & Comments**



IEC 417, No. 5031  
Direct current - Vdc may be used on rating labels.



IEC 417, No. 5032  
Alternating current - For rating labels, the symbol is typically replaced by V and Hz as in 230V, 50Hz. DO NOT USE Vac.



IEC 417, No. 5033  
Both direct and alternating current.



IEC 617-2 No. 02-02-06  
Three-phase alternating current.



IEC 417, No. 5017  
Earth (ground) terminal - Primarily used for functional earth terminals which are generally associated with test and measurement circuits. These terminals are not for safety earthing purposes but provide an earth reference point.



IEC 417, No. 5019  
Protective conductor terminal - This symbol is specifically reserved for the protective conductor terminal and no other. It is placed at the equipment earthing point and is mandatory for all grounded equipment .



IEC 417, No. 5020  
Frame or chassis terminal - Used for points other than protective conductor and functional earth terminals where there is a connection to accessible conductive terminals to advise the user of a chassis connection.



IEC 417, No. 5021

Equipotentiality - Used in applications where it is important to indicate to the operator that two or more accessible functional earth terminals or points are equipotential - More for functional rather than for safety purposes.



IEC 417, No. 5007

On (Supply) - Note that this symbol is a bar, normally applied in the vertical orientation. It is not the number 1.



IEC 417, No. 5008

Off (Supply) - Note that this symbol is a true circle. It is not the number 0 or the letter O.



IEC 417, No. 5172

Equipment protected by double insulation or reinforced insulation (equivalent to Class II if IEC 60536).



ISO 3864, No. B.3.6 Background colour - yellow; symbol and outline – black

Caution, risk of electric shock - Generally used only for voltages in excess of 1000 V. It is permissible to use it to indicate lower voltages if an explanation is provided in the manual. Colour requirements do not apply to markings on equipment if the symbol is moulded or engraved to a depth or raised height of 0.5 mm, or that the symbol and outline are contrasting in colour with the background.



IEC 417, No. 5041 Background colour - yellow; symbol and outline – black

Caution, hot surface - Colour requirements do not apply to markings on equipment if the symbol is moulded or engraved to a depth or raised height of 0.5 mm, or that the symbol and outline are contrasting in colour with the background.



ISO 3864, No. B.3.1 Background colour - yellow; symbol and outline – black

Caution - (refer to accompanying documents) used to direct the user to the instruction manual where it is necessary to follow certain specified instructions where safety is involved. Colour requirements do not apply to markings on equipment if the symbol is moulded or engraved to a depth or raised height of 0.5 mm, or that the symbol and outline are contrasting in colour with the background.



IEC 417, No. 5268-a  
In-position of bistable push control.



IEC 417, No. 5269-a  
Out-position of bistable push control .

# Appendix C - 35V Output Module

## AT35V Output Module

When the Adjustable Mode is enabled for this module, the outputs will provide an adjustable output from 5 volts to 35 volts. The pulse width can be set over the standard range of the unit, but the 35 volt output will self limit to approximately 4  $\mu$ s with some droop. There is no change to TTL output mode functionality with this module.

To maintain the highest possible rise time, care must be taken with cabling and termination. Low capacitance cable and 50 ohm termination will provide the fastest rise times without overshoot. Faster rise times can be achieved by increasing the termination resistance, but some overshoot is likely to occur. While the 35 volt output provides a fast, controlled rising edge, the pulse width and falling edge are not tightly controlled. Also, when using the 35V mode, the option will only function if the 'Polarity' is set for "Active High".

### **35V Specifications** - Through a 50 ohm load at 200 Hz

Output	5V - 35V
Setpoint Resolution	10mV
Rise Time	< 30ns
Accuracy	500mV
Max. Frequency (Internal & External)	4000Hz

# Appendix D - Dual Trigger Input

## **DT15 Dual Trigger Module**

This module option allows the “Gate” input to function as a second trigger input. For consistency, the enabling menu for this option is located under the “Trig” menu structure. Once the dual trigger mode is enabled, both the “Gate” and “Trig” inputs can act as trigger inputs.

Adjustments for the “Gate” trigger input are located under the “Gate” menu structure. The voltage threshold level and trigger edge for the “Gate” input can be adjusted from this menu. The “Gate” trigger edge choice is only available when in dual trigger mode.

Once dual trigger functionality is enabled on the unit, each channel can be assigned to either of the trigger inputs. The default trigger source for each channel is the “Trig” input. The trigger source selection is accessed in the secondary channel menus. To access this menu, first press the yellow “Func” button followed by the channel of interest. Continue to press “Func” then the channel button until the menu page with “Ch Gate:” and “TrigSrc:” appears. Use the “Next” button to place the cursor on the “TrigSrc” line and use the up/down arrows to change to the desired trigger source.

## 9520 SCPI Dual Trigger Command Summary

Keyword	Parameter	Std/ New	Comments
:PULSe [0]		Std	Subsystem. Contains commands to control the output pulse generation. Commands without suffix refer to the currently selected logical instrument. See INSTRument subsystem.
:TRIGger		Std	Subsystem. Contains commands to define the Trigger function.
:MODE	DUAL	New	Sets the unit into dual trigger mode.
:PULSe [1/2/n]		Std	Subsystem. Contains commands to control the output pulse generation. Valid suffix range depends on the number of channels (ChA-1, ChB-2, etc.) Commands without suffix refer to the currently selected logical instrument. See INSTRument subsystem.
:CTRIGger	GATE/TRIG	New	Sets which input is assigned to the channel trigger.

## IEEE 488.2 Common Commands

Mnemonic	Command Name	Parameter	Comments
*TTG	Trigger-Trigger Input		Generates a software trigger pulse for the TRIG input only. Operation is the same as receiving an external trigger pulse.
*GTG	Trigger-Gate Input		Generates a software trigger pulse for the GATE input only. Operation is the same as receiving an external trigger pulse.

# Appendix E - Impedance Matching Outputs

## TZ50 Impedance Matching Output Module

This module option allows a user to have a 50 Ohm load on the output while maintaining an output amplitude of at least 4 Volts while in the TTL/CMOS mode. All other functionality of the module is the same as the AT20 modules, including output while using the Adjustable Mode Function of the channels.

### TTL/ADJUSTABLE OUTPUTS


#### TTL/CMOS Mode

Output Level	4.0v typ into 50 Ohms
Rise Time	3ns
Slew Rate	0.5 V/ns
Jitter - Channel to Channel	50ps RMS

#### ADJUSTABLE MODE

Output Resolution	10mV
Current	200mA typical, 400 mA max (short pulses)
Slew Rate	0.1V/ns

\*Note: The TZ50 module has significant overshoot and ringing through high impedance (see figure).

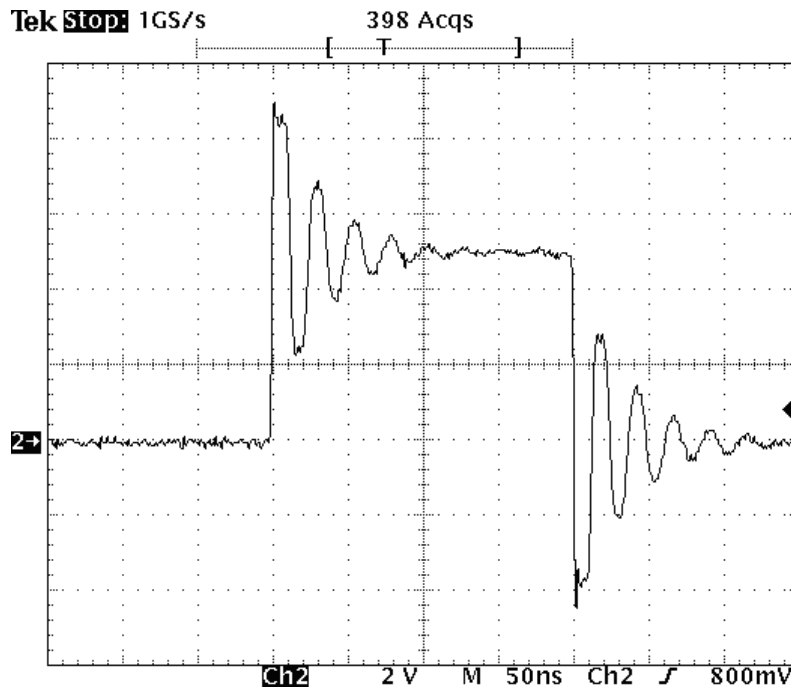

# Appendix F - COM

## Ethernet Interface Overview

An Ethernet interface is optional on the 9520.

The Ethernet module used is a “Digi Connect ME” module supplied by Digi Connectware, Inc. There are several ways to successfully communicate with the pulse generator over Ethernet. The two most popular methods are raw TCP/IP (such as Labview or programming with VISA libraries) and by mapping a PC COM port using the Digi Connectware’s “Realport Drivers”.

Ethernet Communication Notes:

- The Digi Connectware’s “Digi Device Discovery” can be used to determine what IP address was assigned by the local DHCP server (if any).
- “Digi Device Discovery” can also be used to open a web interface to the Ethernet module. Simply double-click on the IP address that is displayed in the Digi Device Discovery utility.

Username: “root”  
Password: “dbps”

- If a mapped COM port is the desired communication method, the Digi Connectware’s “Realport Drivers” setup must be used to install the COM port on the PC. The virtual COM port is then local to the computer it was installed on. Please refer to the Digi Connectware Documentation supplied on the CD, or call Quantum Composers Technical Support.
- The pulse generator’s SCPI parameter “.SYSTEM:COMMunicate:SERial:USB n” is defaulted to 38400 and should not be changed for Ethernet communication, whether or not a mapped COM port is used. The virtual COM port on the PC should be set to 38400 BAUD.
- Echo functionality is not available on the Ethernet port.

# Appendix G - Ethernet Connectivity

## IP Address and Raw TCP/IP Connection

This document describes one of the most popular methods of setting up Ethernet communication for the Quantum Composers pulse generators. The method discussed is Raw TCP/IP communication.


The Ethernet module used in Quantum Composers pulse generators is a “Digi Connect ME” device manufactured by Digi International, Inc. It supports virtually all practical Ethernet communication methods. A set of utilities and documentation by Digi is included on the CD shipped with the pulse generator.

This discussion assumes that the Digi utilities included with your pulse generator and National Instruments VISA (version 3.3 in this procedure, see National Instruments’ website) are installed. The procedures discussed have been prepared using Windows XP service pack 2.

## Determining IP Address

The Digi module has been reset to factory defaults before it left the manufacturing facility. In this mode, it is ready to be assigned an IP address by the local DHCP server. If a crossover cable is being used, the Ethernet device will assume a default IP address.

The Digi utility “Digi Device Discovery” can be used to determine the IP address that is currently assigned to the Ethernet module. Hit “Start, All Programs, Digi Connect, Digi Device Discovery”. When the utility opens, it scans the LAN looking for Digi Ethernet modules. It may take a minute after plugging in or powering the Ethernet module before the LAN negotiates the connection with the Digi module. Hit “Refresh View” in the left column after a minute or so if the utility fails to see the unit when you start it. In some situations it is possible that the Windows Firewall will block the Digi Device Discovery from being able to see the unit. It is advisable to turn the Windows Firewall off while performing these tasks. When the utility sees the Digi device, it will display it in the list (Figure 1).

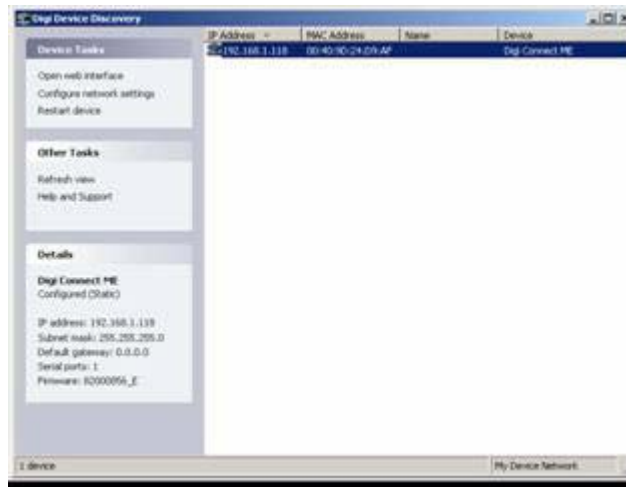

**Figure 1: Digi Device Discovery utility displaying Digi module discovered on the LAN.**

From this point, a web interface can be opened, allowing access to configuration options for the Digi module. If you are required to enter a username and password, they are as follows:

- Username: “root”
- Password: “dbps”

If a static IP address is desired, this change can be made from the web interface. Please note, however, that if the IP address is changed such that it is incompatible with the LAN, all communication including access to the module’s settings (including the IP address!) will no longer be possible over the LAN. If this happens, a crossover cable must be used to access the Digi module’s settings (again using Digi Device Discovery). Temporarily set the pc’s IP address to be compatible with the Digi module’s IP address to get the pc and pulse generator to ‘see’ each other over the crossover cable.

# Appendix H - Increment Mode

## Using the Increment System Mode

The System Increment modes are a pair of special modes, which allow the delay and width of each channel to be incremented at the end of a burst of pulses. Each channel is independent and each may be set with different initial values and different values for the step size for both the delay and the pulse width.

There are two incrementing modes, Increment and DC Increment. In the Increment mode, each start command or external trigger produces a burst of pulses. At the end of the burst the appropriate delays and pulse widths are incremented and the instrument is armed for the next start command. In the DC Increment (Duty Cycle) mode the output is starting as with the normal duty cycle mode. At the end of each cycle the delays and pulse widths are incremented. This continues for the number of cycles defined by the Cycles parameter. The modes are selected from the system mode menu. The step sizes are specified in the channel menus.

## 9520 SCPI Increment Command Summary

Keyword	Parameter	Std/ New	Comments
:PULSe[0]		Std	Subsystem. Contains commands to control the output pulse generation. Commands without suffix refer to the currently selected logical instrument. See INSTRument subsystem.
:MODE	BINCRement /	New	Sets the To mode. Added parameters for Burst Increment and Duty Cycle Increment mode.
:CYCLE	<numeric value>	New	Sets the number of cycles to generate in Duty Cycle Increment mode.
:IRESet	1	New	Resets the width and delay increment parameters on all channels.
:PULSe[1/2/n]		Std	Subsystem. Contains commands to control the output pulse generation. Valid suffix range depends on the number of channels (ChA-1, ChB-2, etc . . .). Command without suffix refers to the currently selected logical instrument. See INSTRuments subsystem.
:IWIDTH	<numeric value>	New	Sets the pulse width increment step size.
:IDELay	<numeric value>	New	Sets the delay increment step size.

**Function-CLR** Pressing the 'FUNCTION' key then 'CLR' initializes the increment parameters and resets the delays and pulse widths to their initial conditions. **This must be pressed after setting all the step parameters, but before generating any pulses.**

#### SPECIFICATIONS

Width Step Size -1.00s to 1.00s

Width Minimum Step 10ns (-10ns)

Width Step Resolution 250ps

Width Incremented Range 10,000s

Delay Step Size -100ms to 100ms

Delay Minimum Step 10ns (-10ns)

Delay Step Resolution 250ps

Delay Incremented Range 10,000s

*Note: Any increment value between -10ns and 10ns will disable the increment function for that parameter.*

Update Rate  $10\mu\text{s} + 30\mu\text{s}$  per active channel  
(1 Ch @ 25 kHz to 8 Ch @ 4 kHz)

# Appendix I - SRS Command Set Compatibility

## Using the SRS Compatible Command Set

Quantum Composers offers SRS Command Set Compatibility as a firmware option to aid customers in transitioning their custom software from the DG535 to the more fully featured QC9520. The majority of these commands function identically to the SRS equivalent instrument capabilities, and several actually extend the functionality and/or the limits of the SRS equivalent capabilities.

		DG535		QC9520	
Command	Example Code	Query Response	Description/Resulting Instrument State	Query Response	Description/Resulting Instrument State
BC {i}		Set or query Pulse/Burst Value between 2 and 32766		Set or query the "On:" Cycles of duty cycle mode (2 - 9,999,999 Pulses). Similar to :SPUL:PCO{?} (see 9520 command reference)	
	CL;BC 3;BC	3	Pulse/Burst = 3	3	on cycles = 3; off cycles = 17
BP {i}		Set or query Pulse/Burst Value between 4 and 32767		Set or query the "Off:" Cycles of duty cycle mode by subtracting the "On:" value from the value received with this command. (4 - 19,999,998 Pulses). Similar to :SPUL:OCO{?} (see 9520 command reference)	
	CL;BP 35;BP	35	Periode/Burst = 35	35	off cycles = 25
CL		Reset to Instrument defaults - #Single Shot Mode, # Set Rate to 10,000 Hz, # 10 pulse per burst, # 20 periods per burst, # Trig Level: +1V DC, # Trig Edge: Rising, # Trig Mode?, # Chan Sync		Set to DG535 Instrument Defaults - difference between DG535 and QC9520 : output amplitudes are set to 2.00V instead of 1.00V	
CS {i}		Set or query Cursor Mode		Unavailable: Always Sets ESR bit 3 - "Wrong Mode for Command"	
DL {i, j, k}		Display Menu Line		Unavailable: Always Sets ESR bit 3 - "Wrong Mode for Command"	
DS {string}		Display a message string of 1 to 20 Characters on the one line display		Display a message string of 1 to 60 characters on the multi-line display	
	DS HELLO WORLD	None	displays "HELLOWORLD"	None	displays "HELLO WORLD"
DT {i, j, t}		Set Delay i linked to j to t seconds ((i or j=1,2,3,4,5,6,7,8,9 => To,A,B,AB,/AB,C,D,CD,/CD)		Set Delay i linked to j to t seconds ((i or j=1,2,3,4,5,6,7,8,9=> To,A,B,C,D,E,F,G,H). Similar to :PULS<i-1>:WIDT{?} in	
	DT 2, 1, 2e-6;DT 2	1, +0.000002000000	Set Delay A to To + 0.000002 seconds	1, +0.000002000000	Set Delay A to To + 0.000002 seconds

		DG535		QC9520	
Command	Example Code	Query Response	Description/Resulting Instrument State	Query Response	Description/Resulting Instrument State
ES {i}		Read and clear entire or specific bit in the Error Status Register		Read and clear entire or specific bit in the Error Status Register	
GT {i} {,j} {,k}		Set 1 to 3 GPIB response terminators		Unavailable: Always Sets ESR bit 3 - "Wrong Mode for Command"	
IC i		Increment or decrement the digit at the current cursor location		Unavailable: Always Sets ESR bit 3 - "Wrong Mode for Command"	
IS {i}		Read and clear entire or specific bit in the Instrument Status Register		Read and clear entire or specific bit in the Instrument Status Register	
MC i		Move cursor left or right one digit		Unavailable: Always Sets ESR bit 3 - "Wrong Mode for Command"	
OA i{,v}		Set or query the output amplitude <v> for channel <i>		Similar to :PULS<i-1>:OUTP:AMPL(?) {v} (see 9520 command reference)	
	OA 3,3,03;OA 3	+3.02	Set Channel B amplitude to +3.02V	+3.02	Set Channel B amplitude to +3.02V
OM i{, j}		Set or query the outputmode <j> of channel <i>		Similar to :PULS<i-1>:OUTP:MOD(?) {mode} (see 9520 command reference)	
	OM 2,3;OM 2	3	Set Channel A output mode to VAR	3	Set Channel A output mode to Adjustable
OO i{,v}		Set or query the output offset <v> for channel <i>		Unavailable: Always sets ESR bit 3 - "Wrong Mode for Command"	
OP i{, j}		Set or query the output polarity <j> for channel <i> (j={0,1}={inverted,normal})		Similar to :PULS<i-1>:POL(?) {polarity} (see 9520 command reference)	
	OP 2,0;OP2	0	Set Channel A output polarity to inverted	0	Set Channel A output to active low
RC i		Recall the instrument settings from location <i> (i=0 is same as CL command => i=0 to 9)		i = 0 to 12 (RC 0 loads the 9520 defaults)	
	RC 0	None	Loads 535 defaults	None	Loads 9520 defaults
	RC 1	None	Load mem location 1	None	Loads mem location 1
SC {i}		Move cursor to column <i>		Unavailable: Always sets ESR bit 3 - "Wrong Mode for Command"	
SM {i}		Set mask for GPIB interrupt service request		Unavailable: Always sets ESR bit 3 - "Wrong Mode for Command"	
SS		Generate a single shot Trigger if Trigger Mode = 2		Similar to *TRG (see 9520 command reference)	
	CL;SS	None	Instrument generates a SS Trigger	None	Instrument generates a SS trigger
ST i		Store instrument settings to location <i>=>1 to 9		i = 0 to 12 for 9520	
	ST 1	None	Stores instrument settings to location 1	None	Stores instrument settings to location 1
TL {v}		Set External Trigger Level to <v> volts		Similar to :SPUL:TRIG:LEV(?) (see 9520 command reference)	
	TL +1.02;TL	+1.02	Set trigger threshold to +1.02V	+1.02	Set trigger threshold to +1.02V
TM {i}		Set Trigger Mode to Int, Ext, SS, or Burst (i=0,1,2,3)		Similar to :SPUL:MOD(?) in combination with :SPUL:TRIG:MOD(?)	
	TM 1;TM	1	Set to Ext Mode	1	Set to Single Shot Mode, Trig Enabled
TR i{,f}		Set Int Trigger Rate (i=0) or Burst Trigger Rate (i=1) to <f>		For better accuracy use :SPUL:PER(?) (see 9520 command reference)	
	TR 1,1e-3;TR 1	0.001		0.001	
TS {i}		Set Trigger Slope to falling or rising edge (i=0,1)		Similar to :SPUL:TRIG:EDG(?) (see 9520 command reference)	
	TS 1;TS	1	Set Trigger to Rising Edge	1	Set Trigger to Rising Edge
TZ i{, j}		Set impedance for Trigger or Outputs (i=0, 1-9) to 50 Ohm or High Z (j=0,1)		Unavailable: Always sets ESR bit 3 - "Wrong Mode for Command"	

# Appendix J - AT45 Option (45 Volt Output)

## Using the AT45 Option

For channels with AT45 output option, the maximum frequency is limited to 1 MHz. The pulse width can be set over the standard range of the unit with both active high and low outputs when set to high impedance mode. In low impedance mode, the pulse width is limited to a maximum of 10s and the active high output is no longer allowed. To maintain the highest possible rise time, care must be taken with cabling and termination. Low capacitance cable and 50 ohm termination will provide the fastest rise times without overshoot. The channel menu structure for the AT45 module changes are described in the table below (changes from standard outputs are in bold italics).

### Channel Menus

Timing Parameters	Pulse Configuration	Pulse Configuration
Channel Enable	Channel Enable	Channel Enable
Sync Source	Polarity: Active High or Active Low	
Pulse Width	<b>Output Type: High Z</b>	<b>Output Type: Low Z</b>
Pulse Delay	<b>Output Level: 3.00 - 45.00 V</b>	<b>Output Level: 3.00 - 45.00 V</b>

Channel Mode	Channel Mode	Channel Mode	Channel Mode	Wait Menu
Channel Enable	Channel Enable	Channel Enable	Channel Enable	
Mode: Normal	Mode Single Shot	Mode: Burst	Mode: Duty Cycle	
			# On Pulses	
		# Pulses in Burst	# Off Pulses	Wait Count